

OCEAN OPTICS UV/VIS SPECTROMETERS INSTRUCTIONS FOR START-UP/SHUT

DOWN—Mac Computers

Updated November 14, 2017

Celeste Morin, x3756, Druckenmiller 243

- 1) Plug the power cable into the wall. The spectrometer fan may turn on at this point. If not, it will turn on in Step #2.
- 2) Plug the USB cable into the spectrometer and back of computer. Fan should be on at this point, if not see your lab instructor.
- 3) Open the Logger Pro software. You should hear a "click" sound when logging in and out of the software. Communication has been established.
- 4) When finished using the spectrometer, unplug the power cable from the wall and unplug the USB cable from the spectrometer. Use the "apple/Q" command to quit out of the Logger Pro software.
- 5) The next set of students must perform Steps 1 through 4.